


# Hydrogen sulphide in liquids and gases

Conform to standard DIN 38405-27

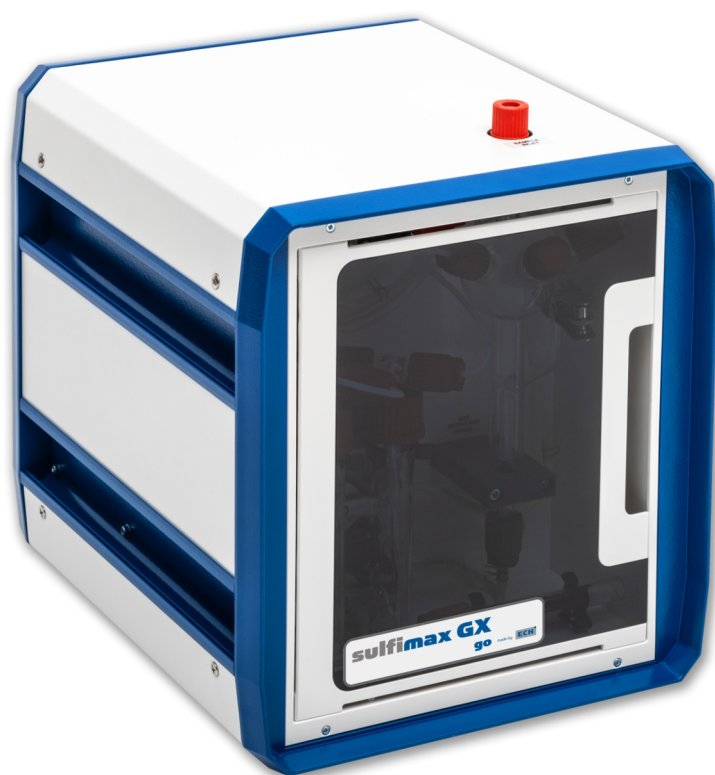

**sulfimax GX**  
**go**

made by **ECH<sup>i</sup>**

# sulfixmax GX

## go

## Hydrogen sulphide in liquids and gases

Conform to standard DIN 38405-27

### Product description

The **Sulfixmax GX Go** determines hydrogen sulphide and volatile sulphides in liquids and gases.

The sample can be dispensed directly by syringe without pretreatment.

By effective gas extraction the  $H_2S$  is completely expelled from the sample.


Interferences due to the sample matrix practically do not occur.

The released  $H_2S$  gas is conducted to the highly sensitive sensor, which detects  $H_2S$  in the range of 0.01 to 10,000 ppm.

A typical measurement takes 5 min, depending on the sample composition.

If the **Sulfixmax GX Go** is extended with the optionally available  $H_2S$  Headspace Module, solid and pasty samples can also be measured.

The Sulfixmax GX Go fulfils the requirements of the standard DIN 38405-27: German standard methods for the examination of water, waste water and sludge - Anions (group D) - Part 27: Determination of sulphide by gas extraction method (D 27).

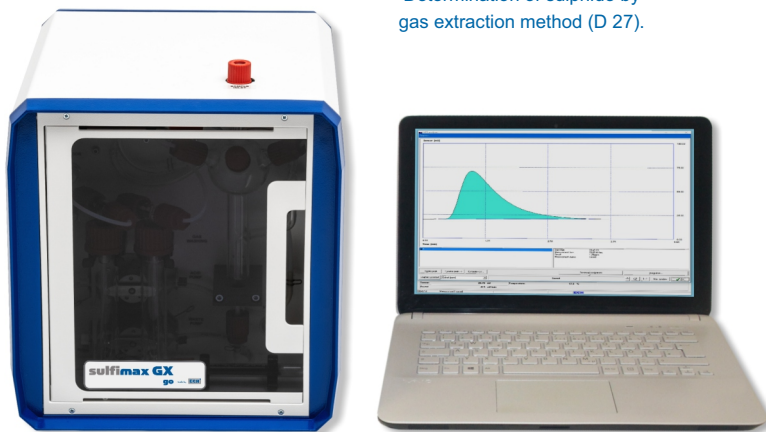


Sulfixmax GX Go - compact version with small footprint

### Applications

- Water, drinking water, surface water
- Municipal wastewater
- Industrial wastewater
- Monitoring of landfill-leachate
- Gas analysis (e. g. LNG, LPG)
- $H_2S$  in hydrocarbon mixtures
- Investigation of technical and pharmaceutical products (e. g. storage stability)
- Quality management

### Advantages



- Analysis of the original sample
- No sample preparation
- Minimized cross sensitivity through the indirect method
- Suitable for on-site use

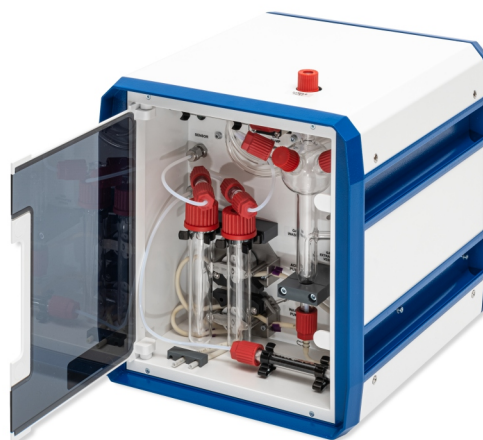

Sulfixmax GX Go - interior view

## Features and Results

- Complete separation of  $\text{H}_2\text{S}$  from the sample by effective gas extraction
- Simple calibration
- Software: simple, clear, intuitive
- Dosing manually by syringe
- Definition of own methods for device control
- Typical measuring time 5 min (depending on sample properties)

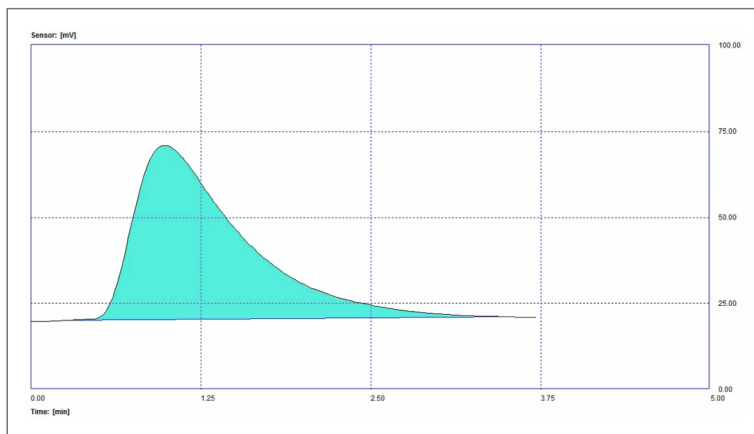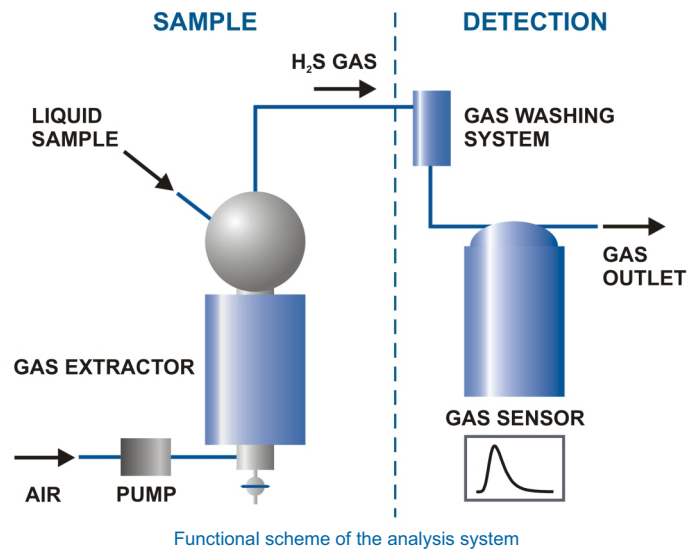

Typical measurement - automatic peak analysis / interpretation

| Result overview |            |                |               |         |
|-----------------|------------|----------------|---------------|---------|
| Mea             | File name  | Sample name    | Sample amount | Result  |
| 1               | 2017110200 | Standard 5 ppm | 1.000 ml      | 5.02 µg |
| 2               | 2017110201 | Standard 5 ppm | 1.000 ml      | 4.81 µg |
| 3               | 2017110203 | Standard 5 ppm | 1.000 ml      | 4.96 µg |
| 4               | 2017110207 | Standard 5 ppm | 1.000 ml      | 4.98 µg |
| 5               | 2017110208 | Standard 5 ppm | 1.000 ml      | 4.93 µg |

|                                 |  |
|---------------------------------|--|
| Evaluation of sub measurements: |  |
| Statistics                      | Arithmetical mean: 4.94 µg<br>Standard deviation: 0.08 µg<br>Rel. standard deviation: 1.58 % |


Value table...

OK Cancel Help

Table of results of a multi measurement

## Technical specifications

|                     |  |
|---------------------|--|
| Measuring range:    | 0.01 ... 10,000 ppm (dependent on sample volume)       |
| Resolution:         | 0.1 µg abs., output signal linear                      |
| Measuring duration: | 2 ... 10 min (dependent on the sample), usually 5 min  |
| Sample volume:      | 0.01 ... 20 mL   |
| Gas flow:           | Up to 50 L/h   |
| Power supply:       | 230 V/50 Hz, 115 V/60 Hz                               |
| Power input:        | 150 W  |
| Dimensions:         | 260 x 310 x 300 mm (W x D x H)                         |
| Weight:             | 8 kg   |
| Data connection:    | RS 232 / USB (with converter)                          |
| Device control:     | PC software (PC not included in the scope of delivery) |

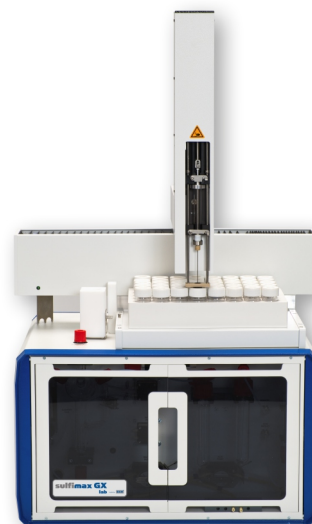

Laboratory version **Sulfimax GX Lab**  
with autosampler for liquids

## H<sub>2</sub>S Headspace Module

### Extension module for solid and viscous samples


The determination of volatile hydrogen sulphide (H<sub>2</sub>S) in solid and viscous samples is easily possible with this module.

It is connected directly to the selective **Sulfimax GX Go** and can be operated by anyone.

The sample is heated isothermally without pretreatment in a sealed headspace vessel and analyzed.

#### Applications:

- Solid samples, e. g. elemental sulfur, sludge, bitumen
- Liquid samples like waste water with sludge particles
- Viscous samples
- Soil samples and waste

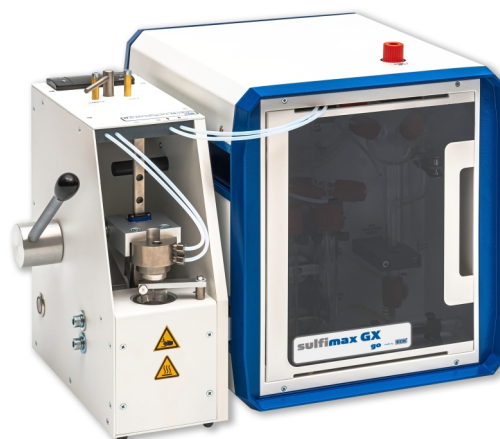

**Sulfimax GX Go** with connected H<sub>2</sub>S Headspace Module

## Book your online demo in the ECH Studio

ECH Scientific have a state of the art laboratory fitted with online presentation capabilities, allowing us to bring product demonstrations live and in full HD, with multiple camera angles and software sharing capabilities enabling us to deliver a full demo experience remotely. Please contact [info@echscientific.com](mailto:info@echscientific.com) to book your session.



ELEKTROCHEMIE HALLE

#### **ECH Elektrochemie Halle GmbH**

*Headquarters in Germany*

Otto-Eißfeldt-Str. 8  
D-06120 Halle (Saale)  
Germany  
+49 (0) 345 279570-0  
[info@ech.de](mailto:info@ech.de)  
[www.ech.de](http://www.ech.de)

#### **ECH Scientific Ltd.**

*Sales and Service Center  
International*

Building 69, Wrest Park, Silsoe  
Bedfordshire, MK45 4HS  
United Kingdom  
+44 (0) 1525 404747  
[info@echscientific.com](mailto:info@echscientific.com)  
[www.echscientific.com](http://www.echscientific.com)

#### **ECH America Inc.**

*Sales and Service Center  
America*

3430 S Sam Houston Pkwy E#400  
Houston TX 77047  
USA  
+1 (713) 805-7525  
[info@echamerica.com](mailto:info@echamerica.com)  
[www.echamerica.com](http://www.echamerica.com)

## the ECH advantage

in-lab | mobile | on-line | process